

BLF Series

Basic series in small scale production, education and research

BLF series is excellent starter and most cost effective equipment for beginners who are new to fermentation in research, education and industrial application. This series requires a laptop / desktop PC for control and operation.

The operation is powered by SOLDAS-BIO software. The graphical operation screen is self explanatory and simple to operate. The BLF series integrated software control measurement and multiple parameter such as temperature, agitation, pumps, pH, DO, etc.

Available in bacterial and cell culture versions, BLF series is supplied as a package with everything necessary to take the next step in capacity and control. For example, the basic bacterial package gives you pH control, feeding, sampling and software for data-logging straight out of the box.




The control system provides all the necessary parameters for a wide range of applications. Common handling issues with standard bench-scale bioreactors have been addressed with interchangeable vessel from 1.5L to 14L, different type of vessel (single wall, double wall, etc) which are removable and autoclavable. The BLF series also offer several ready to use packages and customization request due to price demand.

Ideal For

- Microbial culture growth
(bacteria, yeast and fungi)
- Cell culture growth (Optional Package)
(animal, insect and plant cell)
- Transition from shake | tissue culture flask
- Small scale protein expression
- Education and research

**The images is for Illustration purposes only*

OVERVIEW

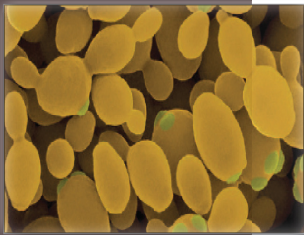

BLF Series

Basic series offer several features to bring flexibility to the user

Key of Features

Base unit:


- Compact, autoclavable laboratory fermentor | bioreactor system
- Multiple choices of heating system (heating block, jacketed vessel, etc)
- Interchangeable single wall vessels with 1.5L, 3L, 5L, 7L, 10L and 14L working volume
- Multiple choice of vessels type (single wall, double wall, etc)
- Interchangeable impeller provide flexibility and high yields for different cell types
- Removable and autoclavable vessel
- Easy handling

Control:

- The system is powered by laptop / desktop pc
- All the measurements and digital controls are controlled by SOLDAS-BIO software
- Adjustable P.I.D values for temperature and pH
- Capability to engage different fermentation mode such batch, fed-batch, and continuous
- Carry out calibration for pumps via SOLDAS-BIO software
- Real time trend data recording and trend display via SOLDAS-BIO software
- Inclusive standard configuration for the system

Standard measurement and control parameters:

- Agitation speed
- Temperature
- pH
- pO₂
- Antifoam / Level
- Feed
- Oxygen supplementation (optional)
- Gas mixing (optional)
- Substrate (optional)
- CO₂ (optional)

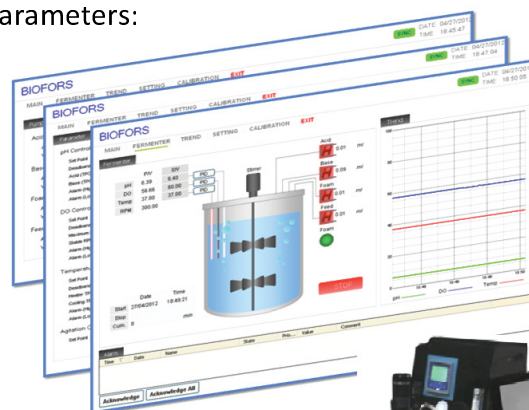

SPECIFICATIONS

Lab Fermenter [BLF-Series]

Culture Vessel

Vessel configuration	1.5L	3L	5L	7L	10L	14L
Total volume	1.5L	3L	5L	7L	10L	14L
Working volume	0.45-1L	0.9-2.1L	1.5-3.5L	2.1-4.9L	3-7L	4.2-9.8L
Material	Borosilicate glass (vessel) SS 316L (top plate)					
Headplate ports	pH DO Anti-foam/ Level Temperature 3/4 way addition Sampling Inoculation Air Inlet Cooler Agitator Exhaust					
Accessories	Exhaust condenser Sparger Cold finger Heating block					
Design	Single-wall or jacketed glass vessel with stainless steel head					

Agitation System

Drive motor	High torque magnetic DC motor, 60 W
Agitation speed	100 ~ 1250 rpm
Driving type	Top driven, Mechanical seal, Direct Coupling
Drive coupling	Direct coupling, easily uncoupled for removal of vessel
Impeller	2x6 blades Rushton turbine (bacterial) / 1x3 blades segment (cell)
Baffle	Vertical three / four plates (removable)

Aeration System

Standard	Air aeration
Microbial (optional)	Air aeration with O ₂ supplementation
Cell (optional)	4-gas mixing system (Air, O ₂ , N ₂ and CO ₂)
Outlet design	Hose tube OD 10mm
Flowmeter	Air callibrated 5 lpm with needle valve and regulator
Inlet Filter	0.2 µm (interchangeable and autoclavable filter)
Sparger / Material	Ring sparger (microbial) µ-sparger (cell culture) / SS 316L
Accuracy	± 5% FS

Peristaltic Pump

Pump Head	Watson Marlow 102R
Motor speed	14.4 rpm ~ 20 rpm
Silicone tube size	0.5 ~ 4.8mm
Max. Flow rate	0.3 ml/min (0.5mm) 0.72 ml/min (0.8mm) 3.07 ml/min (1.6mm) 11.3 ml/min (3.2mm) 22.9 ml/min (4.8mm)
Pump / Control	4 Pumps / On/Off controlled

pH Electrode

Electrode	Mettler Toledo or equivalent
Type	Gel filled, autoclavable
Range	2 ~ 12 pH
Size / Material	Ø12 / Glass
Control	P.I.D controlled

SPECIFICATIONS

Lab Fermenter [BLF-Series]

DO Electrode

Electrode	Mettler Toledo or equivalent
Type	Polarographic, autoclavable
Range	0 ~ 100%
Size / Material	Ø12 / SS 316L
Control	P.I.D controlled

Temperature Electrode

Electrode	BIOFORS
Type	PT100, autoclavable
Range	0 ~ 150°C
Size / Material	Ø4 / SS 316L
Control	P.I.D controlled

Anti Foam Electrode

Electrode	BIOFORS
Type	Conductivity, autoclavable
Holder / Material	Ø6 / Teflon
Controls	Foam control

Reagent Bottle

Reagent	Acid Base Anti-foam Feed bottle
Bottle size	4 x 250 mL
Accessories	0.2 µm PTFE filter (25mm), silicone Tubing

Exhaust Condenser

Type	Double tube water cooled vertical condenser
Material	SS 316L, autoclavable
Accessories	0.2 µm PTFE filter (62mm), silicone Tubing

Requirement

Power supply	120 VAC or 230 VAC 50Hz 10A
Space dimension	850x600x550 (mm) ~ 1000x800x600 (mm); [1.5L ~ 14L]
Water	Controlled @ 2 barg hose connector OD 10mm
Gasses	Controlled @ 1.5 barg , particle and oil free hose connector OD 10mm
Drain	Gravity drain with zero backpressure required hose connector OD 10mm
Laptop (optional)	Windows 7

Malaysia Office:

SOLUTION BIOFORCE SDN. BHD.
 No.3, Jalan TPK 2/4, Taman Perindustrian Kinrara,
 47100 Puchong, Selangor Darul Ehsan, Malaysia
 Tel: +603-80758000 Fax: +603-80755784
 Email: sales@bioforce.com.my
 Web: www.bioforce.com.my

Korea Office:

BIOFORCE CO., LTD.
 Mecha-Zone Na-B126,
 #2979, Hagun-ri, Yangchon-myun,
 Gimpo-si, Gyeonggi-Do,
 415-843 R. O. Korea

